

Special Supplement To Insert Sheet

- When you receive your 100ml sample please use a reduction process to see if your cells can be grown without the presence of FBS. In most cases LiforCell can be used as the only supplement.
- Please use the following example to test LiforCell and FBS combination:
 - If 15% FBS is in use now then the first sample test will use 10% FBS and 5% LiforCell in combination.
 - Then reduce in sample number two to 5% FBS and 10% LiforCell in combination.
 - Then the final sample will be 15% LiforCell and no FBS.

The test should show that LiforCell with your cell will probably require no FBS or less than 5% FBS. To reduce to less than 5% FBS testing should be done at a reduction of 1% FBS and addition of 1% LiforCell until the desired outcome is achieved.

***Lifeblood* Medical, Inc.**

400 Okerson Road, Suite 2B, Freehold, NJ 07728

Phone: 732.431.5833 / Toll Free: 866.389.3335

info@lifebloodmedical.com / www.lifebloodmedical.com